

HDMI FULL HD DISPLAY (13.3") FOR MIKROCAM PRO HDMI AND MINI HDMI MANUAL Art. No. 5914120

A The connecting plate for connection to the microscope camera* is located on the folding support of the display

Accessories:
 2x Mounting screws
 1x Allen (hey) key
 1x HDMI cable
 1x Y-cable hollow plug/USB-C

B The connection plate to the display is located on the side of the microscope camera*

C Attach the display to the microscope camera* and the microscope head*

Fasten to the outside of the folding support using the two mounting screws

D Connect HDMI display via HDMI cable and establish power supply (use camera power supply* (connection 3))

Left-sided interfaces

- 1a Not used
- 2x USB-C
- 2a HDMI

Right-sided interfaces (e.g. MikroCam MINI FULL HD HDMI):

- 2b HDMI
- 1b DC 12V

Connections shown: 2a (HDMI cable), 2b (Y-cable hollow plug/USB-C), 3 (Y-cable hollow plug/USB-C).

E Use the control buttons to adjust the display optimally to your needs

Right-sided interfaces and control buttons

- 4 Menu navigation
- 5 Long: Display ON/OFF
Short: Menu (brightness, contrast)
Not required (USB)

Manual/Accessories/Notes:
www.bresser.de/P5914120

*Not included in the scope of delivery

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail: service@bresseruk.com · Telephone: +44 1342 837098
 BRESSER UK Ltd. · 3G Eden House · Enterprise Way · Edenbridge · Kent TN8 6HF · Great Britain