Digitales Handmikroskop
Digital Handheld Microscope

- DE** BEDIENUNGSANLEITUNG
- EN** INSTRUCTION MANUAL
- FR** MODE D'EMPLOI
- NL** HANDLEIDING
- ES** MANUAL DE INSTRUCCIONES
- IT** MANUALE DI ISTRUZIONI

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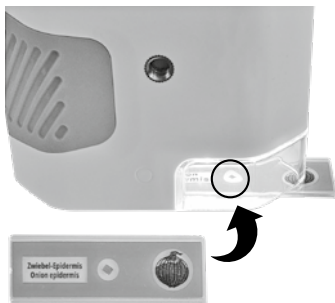
When removing the MicroSD card, gently press the card and wait for the card to pop out. Speed class: C10 or higher. After inserting the card into the device, the device will set this MicroSD card as the preset storage device, the data in the original card may not be readable anymore. Please make a backup of the data first. If the card is not compatible with device, please use the format option in the device, to format the card correctly or change to another card.

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TABLE OF CONTENTS

1. About this manual	10
2. General safety instructions	10
3. Scope of delivery	11
4. Features	11
5. Parts overview	13
6. Charging the device	13
7. Memory Card installation	13
8. Using the device	14
9. Technical data	14
10. Cleaning and Maintenance	15
11. Disposal	15
12. EC Declaration of Conformity	15
13. Warranty & Service	15

WARNING! Not suitable for children under 3 years of age. Small parts, choking hazard. Contains functional sharp edges and points!

WARNING! Only suitable for children aged 8 years and above. Instructions for parents or other responsible persons are included and must be observed. Keep the packaging as it contains important information.

1. ABOUT THIS MANUAL

This instruction manual is to be considered part of the device. Before using the device, please read the safety instructions and the user manual carefully. Keep this user manual for future reference. If the device is sold or passed on, the instruction manual must be

provided to the new owner/user of the product.

2. GENERAL SAFETY INSTRUCTIONS

- **WARNING!** Not for children under three years. Choking hazard – small parts. Contains functional sharp edges and points!
- **ATTENTION:** Only suitable for children of at least 8 years of age. Instructions for parents or other responsible persons are enclosed and must be followed. Keep packaging as it contains important information.
- **RISK OF CHOKING!** Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard.

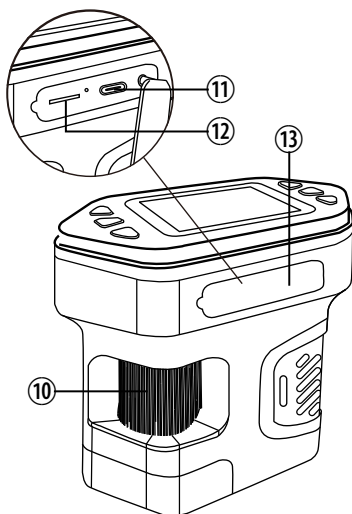
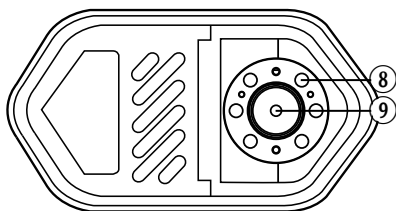
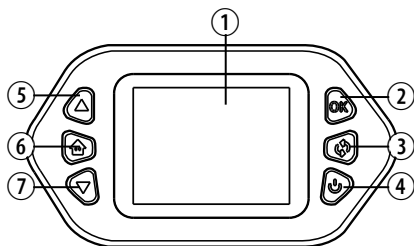
- **RISK OF ELECTRIC SHOCK!** This device contains electronic components that operate via a power source (batteries). Children should only use the device under adult supervision. Only use the device as described in the manual; otherwise, you run the risk of an electric shock.
- **RISK OF CHEMICAL BURN!** Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.
- **RISK OF FIRE/EXPLOSION!** Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands or with different capacities. Remove the batteries from the unit if it has not been used for a long time.
- The manufacturer is not liable for damage related to improperly installed batteries!
- Do not use rechargeable batteries (accumulators).
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Check this device, cables and connections for damage before use.
- Do not immerse the unit in water.
- Do not subject the unit to excessive force, shock, dust, extreme temperature or high humidity, which may result in malfunction, shorter electronic life span, damaged batteries and distorted parts.
- Place your device so that it can be disconnected from the power supply at any time. The power outlet should always be near your appliance and should be easily accessible, as the plug of the power cord serves as a disconnect device from the mains supply.
- To disconnect the unit from the mains, always pull the mains plug and never pull the cable!

3. SCOPE OF DELIVERY

Specimens, carrying strap, USB-C charging cable

4. FEATURES

- Photo, Video, Playback
- Screen size: 2" (5,08 cm) Magnification: Optical: 7x-31x (Digital: 21x-94x)
- Image resolution: 3840x2160P, 1920X1080P
- Video resolution: 1280x720P
- Image format: JPG / Video format: AVI
- Transmission interface: USB Type-C
- Dimensions: 100x95x51 mm / Weight: 157 g

5. PARTS OVERVIEW

1. LCD Screen
2. OK button (Photo/video) 
When the device is in photo mode, press the OK button to take pictures, while the device is in video mode, press the OK button to record video.
3. Mode change 
Press the button several times to choose between different brightness levels or to switch off the lamp.
4. Power button 
Long press the Power button to turn the device on or off.
5. UP button 
In menu options, press the UP button to navigate menu settings upwards. In Photo or video mode, press the UP button to zoom in.
6. Menu button 
Press to open the menu bar with photo / video / playback and system settings. Press the UP/DOWN button to select the desired mode and then press the OK button to confirm.
7. DOWN button 
In menu options, press the DOWN button to navigate menu settings downwards. In Photo or video mode, press the DOWN button to zoom out.
8. LED Lights
9. Microscope lens
10. Focus wheel

11. Charging port
Please use a 5V-1A Power supply to charge the device.
12. MicroSD-card slot
13. Cover charging port

6. CHARGING THE DEVICE

Please do not use the USB port of the computer to directly charge the device. It will automatically switch to the PC CAM mode. To realize the OTG function, you must first turn on the computer, then plug in the USB, and select the desired function. Under the PCCAM mode, you can switch the lens and switch the LED light.

The device has a built-in lithium polymer battery, which allows the device to be used continuously for 2.5 hours after a full charge. The charging time is about 1 hour. The red indicator light is on when charging and goes off when fully charged.

7. MEMORY CARD INSTALLATION

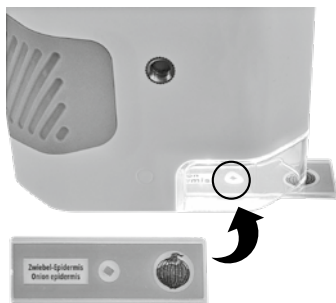
Behind the cover (13) is the MicroSD-card slot. Insert a MicroSD (TF) card into the slot in the correct orientation until the card is fixed in the slot.

When removing the MicroSD card, gently press the card and wait for the card to pop out. Speed class: C10 or higher. After inserting the card into the device, the device will set this MicroSD card as the preset storage device, the data in the original card may not be readable anymore. Please

make a backup of the data first. If the card is not compatible with device, please use the format option in the device, to format the card correctly or change to another card.

8. USING THE DEVICE

Take a specimen slide and put it right under the microscope lens (9). Move around until the specimen is visible in the LCD Display. Use the focus wheel to bring the object into focus. To take a photo, press the OK button.

9. TECHNICAL DATA

Display	2.0-inch color screen
Magnification	1000+
Photo resolution	3840*2160P / 1920*1080P
Video resolution	1280*720P
Focusing roller	Manual focus (0-46mm)
Image format	JPG
Video format	AVI
Dynamic frame rate	20-30 fps
Functions	Photo, Video, Playback, Settings
Transmission method	USB Type-C
Size	100x80.2x50.6 mm
Net weight	156g

10. CLEANING AND MAINTENANCE

Before cleaning the device, disconnect it from the power supply!

Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.

Protect the device from dust and moisture.

11. DISPOSAL

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per Directive 2012/19/EU of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

12. EC DECLARATION OF CONFORMITY

Hereby, Bresser GmbH declares that the equipment type with article number 16443 is in compliance with Directive: 2014/53/EU. The full text of the EU declaration of conformity is available at the following internet address: http://www.bresser.de/download/16443/CE/16443_CE.pdf

13. WARRANTY & SERVICE

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at www.bresser.de/warranty_terms.

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BRESSER Benelux B.V.
Bresser UK Ltd.