MAGUS DIGITAL CAMERA
CDF10 | CDF30 | CDF50 | CDF70

MAGUS

www.magusmicro.com

Levenhuk Inc. (USA):
928 E 124th Ave. Ste D, Tampa, FL 33612, USA,
+1-813-468-3001,
contact_us@levenhuk.com

Levenhuk Optics s.r.o. (Europe):
V Chotejně 700/7, 102 00 Prague 102, Czech Republic,
+420 737-004-919,
sales-info@levenhuk.cz

MAGUS® is a registered trademark of Levenhuk, Inc.

© 2006–2024 Levenhuk, Inc. All rights reserved.

Index

EN	User Manual	4
BG	Ръководство за потребителя	7
CZ	Návod k použití	10
DE	Bedienungsanleitung	13
ES	Guía del usuario	16
HU	Használati útmutató	19
IT	Guida all'utilizzo	22
PL	Instrukcja obsługi	25
PT	Manual do usuário	28
RU	Инструкция по эксплуатации	31
TR	Kullanım kılavuzu	34

EN MAGUS CDF MICROSCOPE DIGITAL CAMERAS

Never look at the sun, a bright light source, or laser beam through the camera. It can cause permanent damage to your eyes.

MAGUS digital cameras are designed to be used with all types of microscopes: biological, fluorescent, metallurgical, polarizing, and stereoscopic. The image of the observed object is displayed on a PC screen. The camera software helps you view, save, and edit captured images and video.

CAMERA INSTALLATION

1. Install the MAGUSView software on your computer from the USB flash drive (included).
2. Connect the camera to your computer using the USB cable.
3. Run the MAGUSView software. The camera will be added in the list of connected devices.
4. Select the camera to start operation.

SPECIAL FEATURES

- The CDF cameras are suitable for darkfield and brightfield microscopy techniques. They are highly light-sensitive and low-noise.
- There are 4 cameras in the CDF series with resolutions ranging from 2 to 20 MP. The higher the magnification of the objective is, the lower the number of pixels that will be required. The extra pixels will have no effect on the image resolution. When working with objectives with magnification up to 40x, however, go for a higher resolution camera. Such a camera captures the details better.
- The frame rate ranges from 20 to 125 fps at the camera's maximum resolution. For studying moving objects, choose a camera with a high frame rate.
- The sensor size is an important parameter for professional shooting. A larger sensor size offers several advantages: cleaner image, more hues, larger field of view. The CDF series offers cameras with sensor sizes from 1/2" to 1".
- The larger the pixel, the more light it collects, which means less image noise and higher light sensitivity of the camera. The CDF series features cameras with pixel sizes ranging from 2.4x2.4 μm to 5.8x5.8 μm .
- The camera outputs the image on the screen in real time. The camera's frame rate is high-speed, allowing you to smoothly adjust the focus when observing with the recommended objective.
- Detailed images with realistic colors are produced when the camera is paired with the right objective.
- The camera is mounted into the vertical tube or a side camera port of the microscope with the C-mount adapter included in the microscope kit. The C-mount adapter (not included) and other adapter rings that match the tube diameter (not included) are used for mounting the camera into the eyepiece. The magnification of the adapter is matched to the size of the camera sensor. The main requirement for adapter optics is to provide a maximum field of view without distortion.
- With the camera software, you can take photos, records video, and edit them. Once calibrated with a calibration slide, the software allows for linear and angular measurements.
- The USB3.0 transfer speed is over 10 times faster than USB2.0. The image is displayed online on the PC monitor. Such speed combined with 30 fps and a higher frame rate transmits the image of moving objects without jerks and delays, and the focus adjustment is smooth even when working with a 100x objective.

THE KIT INCLUDES

MAGUS digital camera, USB cable, USB flash drive with drivers and software, user manual, and warranty card.

SYSTEM REQUIREMENTS

Windows 8/10/11 (32 and 64 bit), Mac OS X, Linux, up to 2.8GHz Intel Core 2 or higher, minimum 2GB RAM, USB 3.0 port.

SPECIFICATIONS

	MAGUS CDF10	MAGUS CDF30	MAGUS CDF50	MAGUS CBF70
Maximum resolution, px	1920x1080	3840x2160	1920x1080	5440x3648
Megapixels	2	8.3	2.1	20
Sensor	1/2" (7.2x4.05mm) SONY Exmor CMOS	1/1.2" (11.14x6.26mm) SONY Exmor CMOS	1/1.2" (11.14x6.26mm) SONY Exmor CMOS	1" (13.06x8.76mm) SONY Exmor CMOS
Pixel size, µm	3.75x3.75	2.9x2.9	5.8x5.8	2.4 x2.4
Light sensitivity	2350mv with 1/30s	5970 mv/lux/s	8935mv with 1/30s	462mv with 1/30s
Exposure	0.1ms–15s	0.02ms–15s	0.014ms–15s	0.1ms–15s
Video recording	+	+	+	+
Frame rate, fps at resolution	125@1920x1080	45@3840x2160 70@1920x1080	96@1920x1080	20@5440x3648 48@2736x1824 58@1824x1216
Image format	*.jpg, *.bmp, *.png, *.tif	*.jpg, *.bmp, *.png, *.tif	*.jpg, *.bmp, *.png, *.tif	*.jpg, *.bmp, *.png, *.tif
Video format	*.wmv, *.avi	*.wmv, *.avi	*.wmv, *.avi	*.wmv, *.avi
Shutter type	ERS	ERS	ERS	ERS
Output	USB3.0, 5Gb/s	USB3.0, 5Gb/s	USB3.0, 5Gb/s	USB3.0, 5Gb/s
Body	metal	metal	metal	metal
Operating temperature range, °C	–10... 50	–10... 50	–10... 50	–10... 50

The manufacturer reserves the right to make changes to the product range and specifications without prior notice.

CARE AND MAINTENANCE

- Do not disassemble the camera on your own.
- Keep the camera away from moisture; do not use it in rainy weather.
- Keep the camera away from shock or excessive pressure.
- Do not overtighten the locking screws.
- Keep the camera away from hazardous environment, home and car heaters, incandescent lamps or open fire.
- When cleaning the lenses, first blow any dust or debris off the surface or wipe them off with a soft brush. Then wipe the lens with a soft cleaning cloth slightly damp with spirits or ether.
- Seek medical advice immediately if a small part or a battery is swallowed.

MAGUS WARRANTY

Magus optic products carry a **5-year warranty** against defects in materials and workmanship. All Magus accessories are warranted to be free of defects in materials and workmanship for **six months** from the purchase date. The warranty entitles you to the free repair or replacement of the Magus product in any country where a Levenhuk office is located if all the warranty conditions are met.

For further details, please visit: www.magusmicro.com

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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