
- DE Bedienungsanleitung**
- GB Operating Instructions**
- FR Mode d'emploi**
- NL Handleiding**
- IT Istruzioni per l'uso**
- ES Instrucciones de uso**

SERVICE AND WARRANTY:

www.bresser.de/warranty_terms

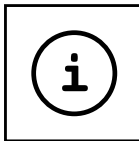MICROSCOPE GUIDE:

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MANUAL DOWNLOAD:

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
MICROSCOPE FAQ:

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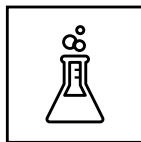


EXPERIMENTS:

www.bresser.de/downloads

DE	Bedienungsanleitung.....	4
GB	Operating Instructions	10
FR	Mode d'emploi.....	16
NL	Handleiding	22
IT	Istruzioni per l'uso	28
ES	Instrucciones de uso	34

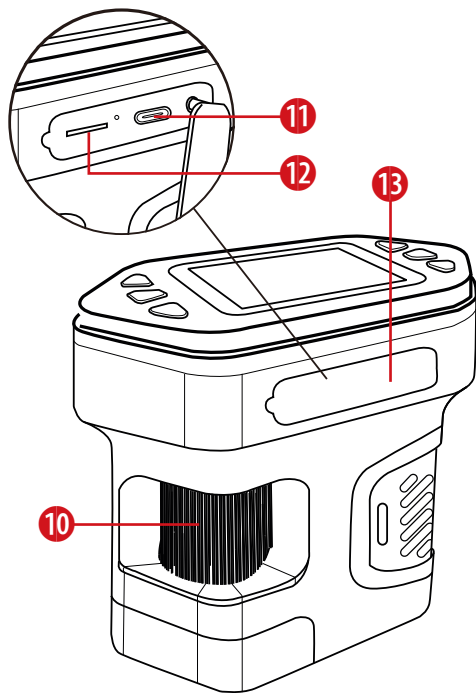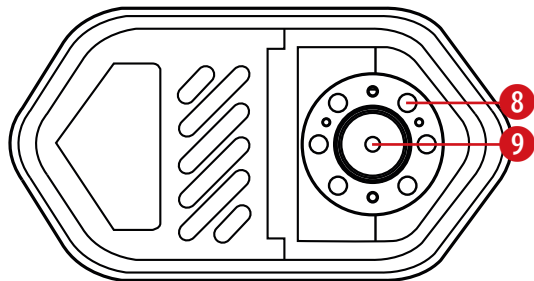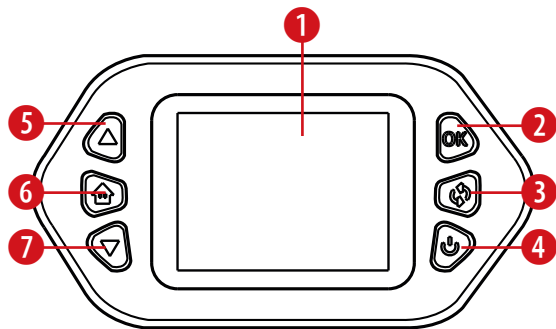DE

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EN

Operating Instructions

WARNING! Not for children under three years. Choking hazard – small parts. Contains functional sharp edges and points!

ATTENTION: Only suitable for children of at least 8 years of age. Instructions for parents or other responsible persons are enclosed and must be followed. Keep packaging as it contains important information.

About this Instruction Manual

Please read the safety instructions and the operating instructions carefully before use. Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided

to the new owner/user of the product.

General Warnings

- **Risk of choking** – Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard. This product contains small parts that could be swallowed by children.
- Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.
- **Risk of electric shock** – Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections

for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

- **Risk of property damage!** – Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose the device to high temperatures. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- Do not interrupt an on-going data connection to a computer before all data has been transferred. This could lead to data loss for which the manufacturer is not liable.
- This toy contains batteries that are non-replaceable.

Notes on cleaning

- Turn off the device
- Clean the lenses only with a soft, lint-free cloth, (e.g. microfibre cloth). To avoid scratching the lenses, use only gentle pressure with the cleaning cloth.
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.

Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per the Directive 2012/19/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German

law, used electronic devices must be collected separately and recycled in an environmentally friendly manner. Empty, old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 6 January 2006 from your local waste-disposal service or environmental authority.

Drain the battery completely before disposal.



EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address:
www.bresser.de/download/8855200/CE/8855200_CE.pdf

Scope of delivery

Digital Hand Microscope, Specimens, Carrying strap, USB-C Charging cable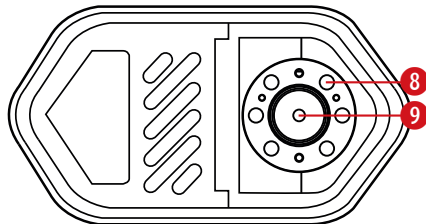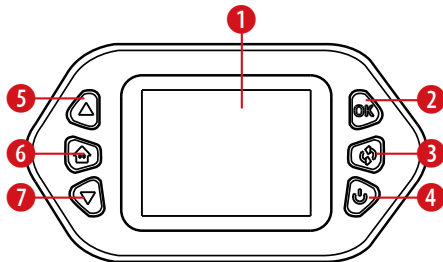

Parts overview

1. LCD Screen

2. OK button (Photo/video)

When the device is in photo mode, press the OK button to take pictures, while the device is in video mode, press the OK button to record video.

3. Mode change

Press the button several times to choose between different brightness levels or to switch off the lamp.

4. Power button

Long press the Power button to turn the device on or off.

5. UP button

In menu options, press the UP button to navigate menu settings upwards. In Photo or video mode, press the UP button to zoom in.

6. Menu button

Press to open the menu bar with photo / video / playback and system settings. Press the UP/DOWN button to select the desired mode and then press the OK button to confirm.

7. DOWN button


In menu options, press the DOWN button to navigate menu settings downwards. In Photo or video mode, press the DOWN button to zoom out.


8. LED Lights

9. Microscope lens

10. Focus wheel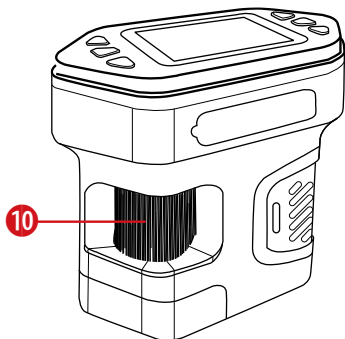

Charging the device

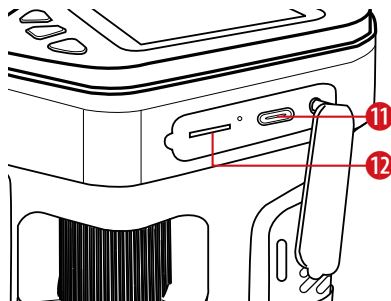10. Charging port: Please use a 5V-1A Power supply to charge the device.

Please do not use the USB port of the computer to directly charge the device. It will automatically switch to the PC CAM mode. To realize the OTG function, you must first turn on the computer, then plug in the USB, and select the desired function. Under the PCCAM mode, you can switch the lens and switch the LED light.

The device has a built-in lithium polymer battery, which allows the device to be used continuously for 2.5 hours after a full charge. The charging time is about 1 hour. The red indicator light is on when charging and goes off when fully charged.


Memory Card installation

11. Behind the cover (13) is the MicroSD-card slot. Insert a MicroSD (TF) card into the slot in the correct orientation until the card is fixed in the slot.

When removing the MicroSD card, gently press the card and wait for the card to pop out. Speed class: C10 or higher. After inserting the card into the device, the device will set this MicroSD card as the preset storage device, the data in the original card may not be readable anymore. Please make a backup of the data first. If the card is not compatible with device, please use the format option in the device, to format the card correctly or change to another card.

Using the device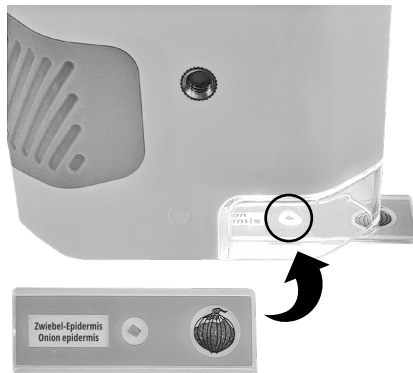

Take a specimen slide and put it right under the microscope lens (9). Move around until the specimen is visible in the LCD Display. Use the focus wheel to bring the object into focus. To take a photo, press the OK button.

Technical Data

Display	2.0-inch color screen
Magnification	1000+
Photo resolution	3840*2160P / 1920*1080P
Video resolution	1920*1080P / 1280*720P
Focusing roller	Manual focus (0-46mm)
Image format	JPG
Video format	AVI
Dynamic frame rate	20-30 fps
Functions	Photo, Video, Playback, Settings
Transmission method	USB Type-C
Size	100x80.2x50.6 mm
Net weight	156g

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Bresser UK Ltd.