7" LCD-Mikroskop
7" LCD Microscope

- DE** **BEDIENUNGSANLEITUNG**
- EN** **INSTRUCTION MANUAL**
- FR** **MODE D'EMPLOI**
- NL** **HANDLEIDING**
- IT** **MANUALE DI ISTRUZIONI**
- ES** **MANUAL DE INSTRUCCIONES**

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TABLE OF CONTENTS

1. About this manual	14
2. General safety instructions	14
3. Scope of delivery	15
4. Features	15
5. Parts Overview	17
6. Component functions and descriptions	18
7. Remote control	18
8. Handheld Microscope	19
9. Photo/Video Storage, Transfer & PC Viewing	20
10. TF Card Installation & Precautions	20
11. Charging Methods	21
12. Charging & Indicator Status	21
13. Battery Usage & Safety Precautions	22
14. Important Precautions	22
15. Product Specifications	22
16. Cleaning and Maintenance	23
17. Disposal	23
18. EC Declaration of Conformity	23
19. Warranty & Service	23

WARNING! Not suitable for children under 3 years of age. Small parts, choking hazard. Contains functional sharp edges and points!

WARNING! Only suitable for children aged 8 years and above. Instructions for parents or other responsible persons are included and must be observed. Keep the packaging as it contains important information.

1. ABOUT THIS MANUAL

This instruction manual is to be considered part of the device. Before using the device, please read the safety instructions and the user manual carefully. Keep this user manual for future reference. If the device is sold or passed on, the instruction manual must be provided to the new owner/user of the product.

2. GENERAL SAFETY INSTRUCTIONS

- **RISK OF CHOKING!** Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard.
- **RISK OF ELECTRIC SHOCK!** This device contains electronic components that operate via a power source (batteries). Children should only use the device under adult supervision. Only use the device as described in the manual; otherwise, you run the risk of an electric shock.
- **RISK OF CHEMICAL BURN!** Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.





- **RISK OF FIRE/EXPLOSION!** Use only the recommended batteries. Do not short-circuit the device or batteries, or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- Use only the recommended batteries. Always replace weak or empty batteries with a new, complete set of batteries at full capacity. Do not use batteries from different brands or with different capacities. Remove the batteries from the unit if it has not been used for a long time.
- The manufacturer is not liable for damage related to improperly installed batteries!
- Do not use rechargeable batteries (accumulators).
- Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Check this device, cables and connections for damage before use.
- Do not immerse the unit in water.
- Do not subject the unit to excessive force, shock, dust, extreme temperature or high humidity, which may result in malfunction, shorter electronic life span, damaged batteries and distorted parts.
- Place your device so that it can be disconnected from the power supply at any time. The power outlet should always be near your appliance and should be easily accessible, as the plug of the power cord serves as a disconnect device from the mains supply.
- To disconnect the unit from the mains, always pull the mains plug and never pull the cable!

3. SCOPE OF DELIVERY

Handheld microscope, base, handle, petri dishes, box with specimens, collection box, tweezers, pipette, USB-C cable

4. FEATURES

1. Optimal student microscope
2. Microscope with XXL screen (7")
3. Photo and video recording possible
4. Microscope can be used as a microscope or monocular
5. Battery operated

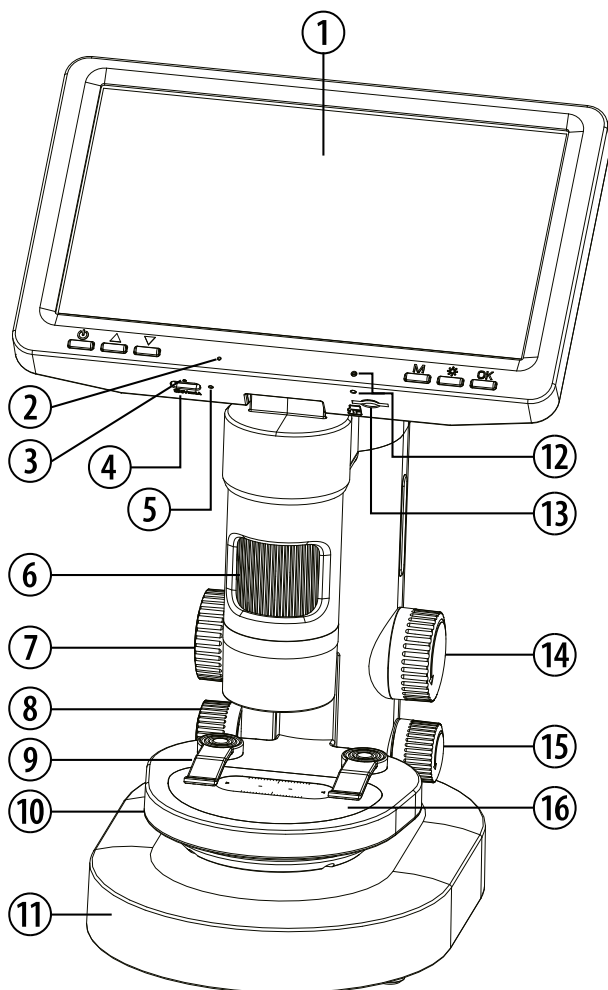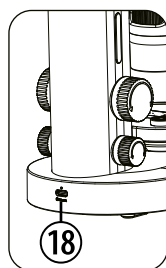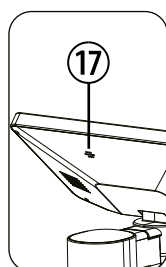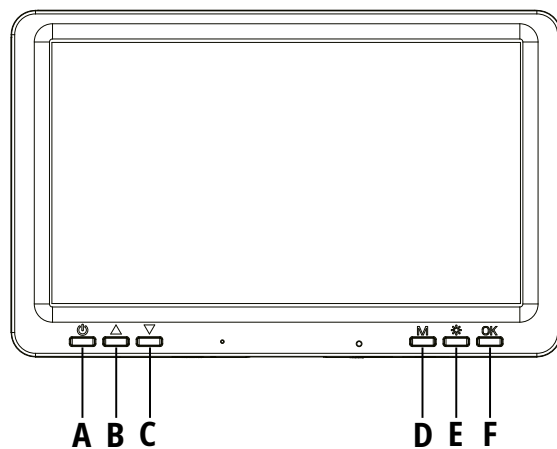5. PARTS OVERVIEW

- A. Power/Delete Button
- B. Up/Magnify Button
- C. Down/Reduce Button
- D. Menu/Return/Flip Button
- E. LED Lens Light Button
- F. Confirm/Filter Button
- 1. Screen
- 2. Microphone hole
- 3. Charging indicator light
- 4. USB Type-C charging port 1
- 5. Reset button
- 6. Focusing wheel
- 7. Stage adjustment wheel
- 8. Observation window switch
- 9. Dust removal cover
- 10. Stage platform
- 11. Base
- 12. Infrared receiver
- 13. TF card slot
- 14. Stage adjustment knob
- 15. Observation window switch
- 16. Observation window
- 17. Speaker
- 18. USB Type-C charging port 2

6. COMPONENT FUNCTIONS AND DESCRIPTIONS

Power/Delete Button

- When powered off: Long press to turn on
- In any interface: Long press to power off
- Image mode: Short press to display “Delete” option

Up/Magnify Button

- In menu: Short press to select “Previous item”
- In observation mode: Short press to zoom in

Down/Reduce Button

- In menu: Short press to select “Next item”
- In observation mode: Short press to zoom out

Menu/Return/Flip Button

- When powered on: Short press to access options (“Photo”, “Video”, “Album”, “Settings”)
- In menu mode: Short press to return
- In observation mode: Long press to vertically flip the interface

LED Lens Light Button


- When powered on: Short press to adjust brightness (cycles through 6 levels + Off)

Confirm/Filter Button

- In menu: Short press to confirm selection
- In capture mode: Short press to take photo/start recording
- In observation mode: Long press (3 seconds) to activate filters

7. REMOTE CONTROL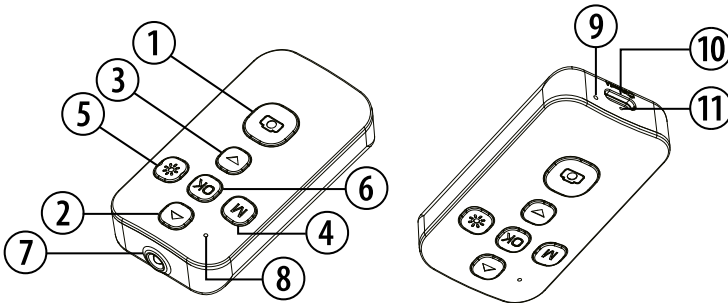

1. Photo/Video button

- Photo mode: Press briefly to take a photo.
- Video mode: Press briefly to record a video.

2. Up/Zoom-in button

- Menu mode: Press briefly to select the previous item in the menu/settings.
- View mode
- Press briefly to zoom in.

3. Decrease/Zoom out button

- Menu mode: Press briefly to select the next item in the menu/settings.
- View mode: Press briefly to zoom out.

4. Menu/Back button

- When switched on: Press briefly to open the options menu: Photo / Video / Album / Settings.
- Menu mode: Press briefly to return to the previous screen.

5. LED light button


- When switched on: Press briefly to adjust the brightness (cycles through the levels: 1, 2, 3, 4, 5, 6, Off).

6. OK/Confirm button


- Menu mode: Press briefly to confirm the selected menu/setting.

7. Infrared transmitter**8. Status indicator****9. Reset hole****10. Type-C charging port****11. Charging indicator****8. HANDHELD MICROSCOPE**

- The microscope main unit can be removed separately

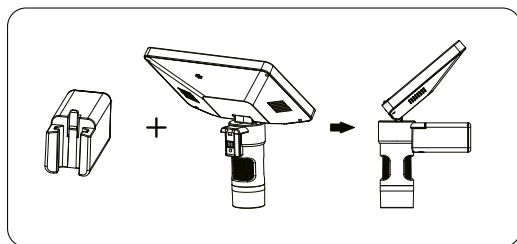
- Attach the microscope main unit to the handle to use as a handheld microscope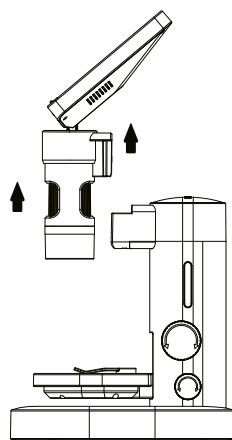

9. PHOTO/VIDEO STORAGE, TRANSFER & PC VIEWING


1. Connect to the computer via a Type-C data cable.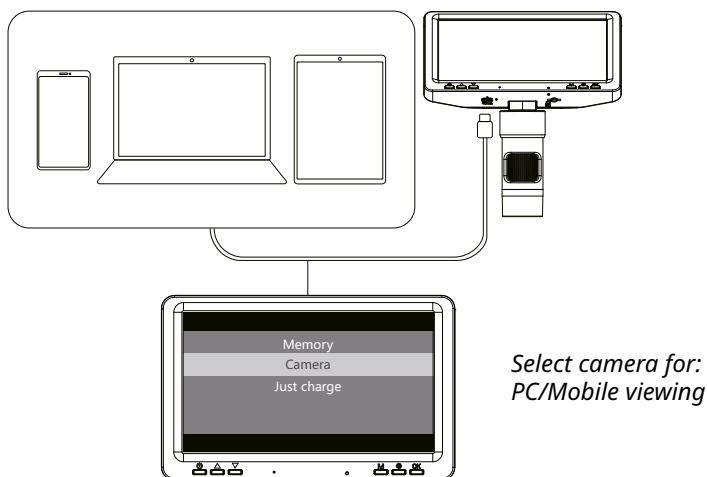

2. Supports data transfer and viewing on select Windows, macOS, and Android systems (iOS and iPadOS not supported)

10. TF CARD INSTALLATION & PRECAUTIONS

1. Insert the TF card into the slot in the correct orientation until it clicks into place. To remove, gently push the card to release it. (Only Class 10 or higher speed TF cards are supported. Never insert/remove the card while the device is operating to prevent damage.)
2. Ensure proper insertion direction. Reverse insertion may damage both the device and TF card slot.
3. When inserted, the device will automatically set the TF card as default storage. Existing data may become inaccessible - back up your files first.
4. Always format the TF card using the device's settings menu before first use. If incompatible, try another TF card.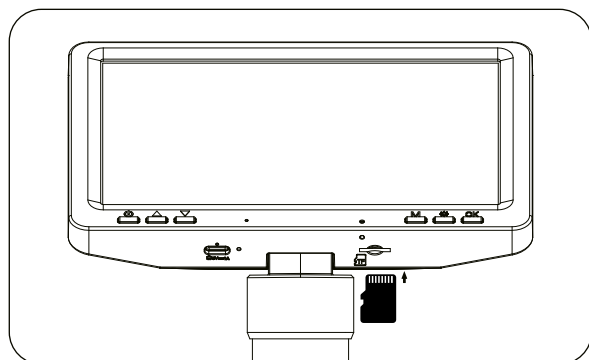


*Installation: Insert it first
and then press in*

*Take it out: Press it first
and then pull it out*


11. CHARGING METHODS

Charging Method 1

The handheld unit and base can be charged separately.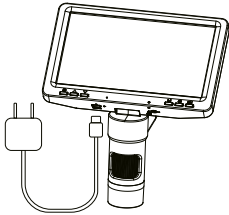

*Type-C Charging Port 1
(Handheld Unit)*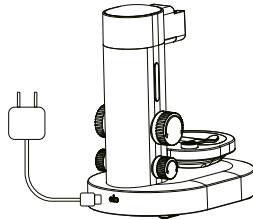


*Type-C Charging Port 2
(Base Unit)*

Charging Method 2

When the base is powered:

- Place the handheld unit on the base stand
- Align power connection points 1 & 2 for system charging

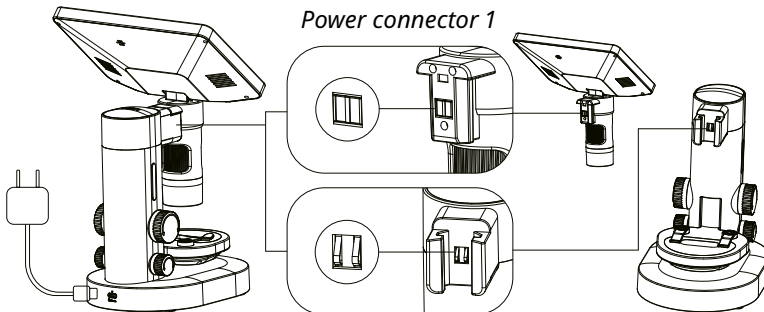Type-C charging port 2

Power connector 2

Note: When the system is charging, You can also charge using the handheld unit's Charging Port 1. Just so you know, the base power connection will automatically disconnect when you use Port 1.

12. CHARGING & INDICATOR STATUS

Charging Instructions:

1. When battery is low, charge immediately using a 5V-1A power adapter. Using the device while charging may reduce battery lifespan.
2. Do not charge via computer USB port while powered on - this automatically activates PC camera mode.
3. Compatible with power banks for charging.

Indicator Lights:

- The red charging indicator light is on, and it will turn off when fully charged.

13. BATTERY USAGE & SAFETY PRECAUTIONS

1. The device contains one built-in 3.7V/1000mAh polymer lithium battery and one 3.7V/3000mAh polymer lithium battery, which are non-removable and non-replaceable. Contact the seller for repairs if damaged.
2. Batteries must be charged under adult supervision.
3. Turn off the device switch when charging the batteries.
4. Do not short-circuit the power terminals.
5. Do not connect to power sources exceeding the recommended quantity or voltage (Recommended adapter specifications: 5V 1A).
6. During use, regularly inspect wires, plugs, casings and other components for damage. Stop use immediately if damage is found and do not resume until fully repaired.
7. When not using the device for extended periods, fully charge the battery before turning off the power switch.

14. IMPORTANT PRECAUTIONS

1. This device is a precision imaging product. Do not knock, impact, squeeze, or drop it.
2. Contains built-in polymer lithium battery. Comply with local battery usage and recycling regulations.
3. The product is not waterproof. Do not immerse in water or store in humid environments.
4. Do not expose the product to direct sunlight for extended periods.
5. Avoid using in high-humidity environments as this may damage the motherboard.
6. After each use, store the product properly. Do not leave it unattended.
7. Do not clean with corrosive cleaners as this may cause damage.
8. If unused for extended periods, charge the device at least once per month.
9. If the device freezes or malfunctions during operation, press the reset button once.

15. PRODUCT SPECIFICATIONS

- Dimensions: 181x170x298.5 cm
- Display: 7-inch color screen
- Magnification: 200x10
- Photo Resolution: 5120x2880 2048x1152
- Video Resolution: 2560x1440 1920x1080
- Focusing Mechanism: Manual focus (0-46 mm)
- Image Format: JPG
- Video Format: mp4
- Frame Rate: 20-30 f/s
- Max Supported Capacity: 128 G

- Functions: Photo Video Microscope Telescope Magnifier
- System Languages: DE, EN, CN, FR, PT, PL, NL, TR
- Interface: Type-C
- Weight: 686 g

16. CLEANING AND MAINTENANCE

Before cleaning the device, disconnect it from the power supply!

Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.

Protect the device from dust and moisture.

17. DISPOSAL

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per Directive 2012/19/EU of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner.

18. EC DECLARATION OF CONFORMITY

Hereby, Bresser GmbH declares that the equipment type with article number 16461 is in compliance with Directive: 2014/53/EU. The full text of the EU declaration of conformity is available at the following internet address:
http://www.bresser.de/download/16461/CE/16461_CE.pdf

19. WARRANTY & SERVICE

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at www.bresser.de/warranty_terms.

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BRESSER Benelux B.V.
Bresser UK Ltd.