FREEK VONK®

×

(B) BRESSER®

DIGITAAL HANDHELD MICROSCOOP
DIGITAL HANDHELD MICROSCOPE
DIGITALES HANDMIKROKSOP
MICROSCOPE PORTATIF NUMÉRIQUE
MICROSCOPIO DIGITAL DE MANO

Art. No. 9820300

- (NL)** Handleiding
- (EN)** Operating Instructions
- (DE)** Bedienungsanleitung
- (FR)** Mode d'emploi
- (ES)** Instrucciones de uso

SERVICE AND WARRANTY:

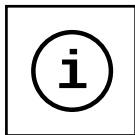
www.bresser.de/warranty_terms

MANUAL DOWNLOAD:

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
MICROSCOPE GUIDE:

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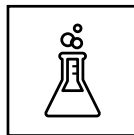EXPERIMENTS:

www.bresser.de/downloads






MICROSCOPE FAQ:

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(NL)	Handleiding.....	4
(EN)	Operating Instructions	10
(DE)	Bedienungsanleitung	16
(FR)	Mode d'emploi	22
(ES)	Instrucciones de uso	28

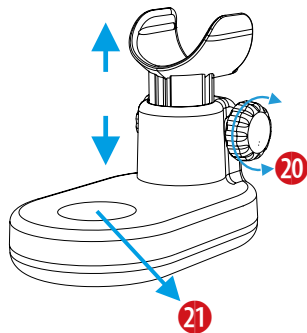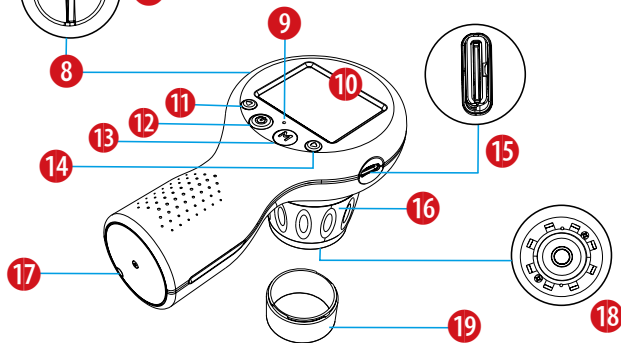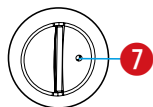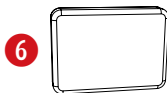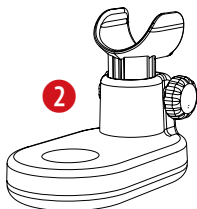Handleiding

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Operating Instructions

WARNING! Not for children under three years. Choking hazard – small parts. Contains functional sharp edges and points!

ATTENTION: Only suitable for children of at least 8 years of age. Instructions for parents or other responsible persons are enclosed and must be followed. Keep packaging as it contains important information.

About this Instruction Manual

Please read the safety instructions and the operating instructions carefully before use. Keep these instructions for renewed use at a later date. When the device is sold or given to someone else, the instruction manual must be provided

to the new owner/user of the product.

General Warnings

- **Risk of choking** – Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard. This product contains small parts that could be swallowed by children.
- Leaking battery acid can lead to chemical burns. Avoid contact of battery acid with skin, eyes and mucous membranes. In the event of contact, rinse the affected region immediately with a plenty of water and seek medical attention.
- **Risk of electric shock** – Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections

for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.

- **Risk of property damage!** – Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose the device to high temperatures. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- Do not interrupt an on-going data connection to a computer before all data has been transferred. This could lead to data loss for which the manufacturer is not liable.

Disposal

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per the Directive 2012/19/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner. Empty, old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices or batteries produced after 6 January 2006 from your local waste-disposal service or environmental authority.

Drain the battery completely before disposal.

EC Declaration of Conformity

Bresser GmbH has issued a "Declaration of Conformity" in accordance with applicable guidelines and corresponding standards. The full text of the EU declaration of conformity is available at the following internet address:
www.bresser.de/download/9820300/CE/9820300_CE.pdf

Scope of delivery

Handheld microscope, base, bag, specimens, carrying strap, USB-C cable

Your microscope is made up of the following parts:

- 1 Handheld Microscope
- 2 Base - For placing leaves, small stones, slides, etc.
- 3 USB Type-C cable (for charging)
- 4 Carrying strap
- 5 Cleaning cloth
- 6 Blank utensil box
- 7 Reset button
- 8 MicroSD/TF card slot
- 9 Charging indicator light - green = full, red = charging
- 10 Display
- 11 LEFT/Digital zoom (short press), Brightness down (long press)
- 12 Confirm (take photos/videos) button (short press), Power on/off (long press)
- 13 MODE (switch between photo/video/playback) button (short press), Menu button (long press)
- 14 RIGHT/Digital zoom (short press), Brightness up (long press)
- 15 USB Type-C charging port

- 16 Focussing wheel
- 17 Carrying strap hole
- 18 LED-Illumination
- 19 Translucent cover - When observing the specimen, you can clearly see the current status. The height of the cover is exactly to the best focussing distance of the microscope), removable
- 20 Table height adjustment
- 21 Microscope table observation area

Location

Before you start, choose an ideal location for using your microscope. It is important that you have enough light available (window, desk lamp).

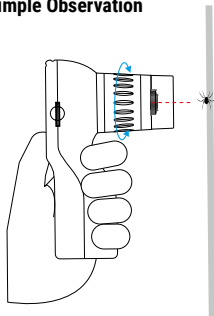
Charging instructions

Insert the USB cable into the microscope USB Type-C port and connect the other end to a USB power supply for charging. Charging time is about 1.5 hours to fully charge. Use time is about 3 hours. The indicator light turns red when charging and green when fully charged.

MicroSD/TF Memory Card

Insert the microSD/TF card into the slot with the back facing upwards.

Simple Observation

Long press the power button (12) to activate the screen. Place the object to be observed at the center of the screen. Short press the LEFT (11) or RIGHT button (14) to adjust the zoom level and use the focussing wheel (16) to adjust picture clarity.

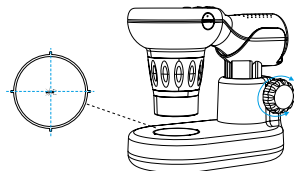
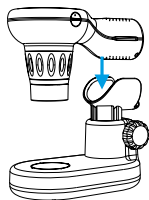
When observing objects, adjust the table height, focussing wheel and zoom level accordingly to obtain a clear image. Avoid excessive zooming in or out.

LED-Illumination

Press and hold the RIGHT (14) button to switch the light (18) on or increase the brightness. Press and hold the LEFT (11) button to reduce the brightness or switch off the light.

Using the base

Snap the handheld microscope handle onto the base (25).

Place the object to be observed in the center of the observation area of the microscope table. Adjust the table height and observe the changes on the screen. When the picture becomes clear, you can take pictures or videos.

Mode switching

Short press the mode button (13) to cycle through the following three Modes: Photo/Video/Playback.

Taking pictures/videos

When in photo or video mode, short press the confirm button (12) to take a photo or start a video.

Hold the microscope steady to avoid blurred images.

Settings mode

Long press the mode button (13) to enter the settings mode. Use the LEFT or RIGHT buttons to navigate the menu. Short press the confirm button (12) to enter menu options.

Playback mode

When playing a video, press and hold the LEFT or RIGHT button to fast-forward or rewind.

Delete files

Select the photo or video in playback mode that you want to delete. Long press the MODE button and choose delete. Press the POWER button to confirm.

Experiments

Use the following web link to find interesting experiments you can try out.

<http://www.bresser.de/downloads>

Troubleshooting

Question	Solution
The microscope won't turn on	Maybe the battery is dead. Please charge.
The picture has a lot of noise and is very dark	Please turn on the light.
System stuck	Please insert a thin needle into the reset hole and press it to reset the system.
The picture is not clear	Please check whether the focus is correct and whether the object being observed is within the focus range.
The picture is too bright or too dark	Please enter the settings menu to adjust the exposure parameters.

TIPS on cleaning

Remove the device from its energy source before cleaning (remove plug from socket).

Clean the exterior of the device with a dry cloth. Do not use cleaning fluids so as to avoid causing damage to electronic components.

Protect the device from dust and moisture. Store the device in the bag supplied or in its original packaging.

Warranty & Service

The regular guarantee period is 2 years and begins on the day of purchase. To benefit from an extended voluntary guarantee period as stated on the gift box, registration on our website is required.

You can consult the full guarantee terms as well as information on extending the guarantee period and details of our services at www.bresser.de/warranty_terms.

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© en onder licentie van Studio Freek,
Bresser UK Ltd.